User and maintenance manual and CE certification

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BENCH FUME HOOD  
Series models AB - serial number:

**prodotto da/produced by**  
IRIDE International srl  
Via S.Giuseppe 5, Azzano Decimo - PN  
Italia

**emesso da/issued by**  
IRIDE International srl  
Via S.Giuseppe 5, Azzano Decimo - PN  
Italia

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### BENCH FUME HOOD GENERAL DESCRIPTION

The bench fume hood is made of a galvanised steel plate case painted with resistant epoxy paint and a 255 Watt induction motor.

The following are the specifications of the filters supplied with it: 9,200 cm<sup>2</sup> nominal filter surface, 20 micron filter nominal capacity; one is a pre-filter located below the motor which does not need maintenance as it is meant solely to protect the motor from any misuse of the other filter which must be regularly cleaned and replaced.

The control panel comprises the following switches:

1 16A suction switch

Suction 550 m<sup>3</sup>/h

Suction is powered by a 220Volt/50Hz voltage.

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*The hood comes with a one-year period warranty which does not cover any damage or misuse caused by the user.*

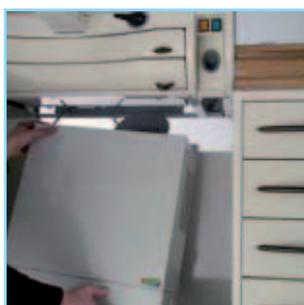
*Use the hood carefully and in a proper way, avoiding any improper or faulty installation breaching this warranty or safety rules in compliance with its manufacturing standards.*

*Follow this manual very carefully, concerning the proper use and maintenance of the bench fume hood.*

*Before using the appliance, any operator must be perfectly trained on its functions and technical specifications.*


*A filter is supplied with the appliance; the package also contains: 1 user and maintenance manual and CE certification.*

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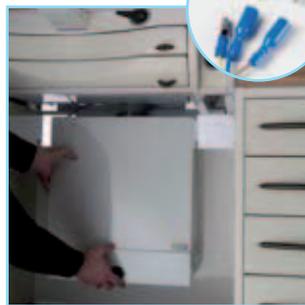*Grasp the hood and insert it (see picture) under the bench, securing it to the bench guide bars located at the back.*

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*Make sure that the hood is completely leant and secured to the back, then connect the electrical system.*

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*Connect the electrical system located under the control panel as shown in the picture.*

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*Finally, insert the ductwork and lock it to the air opening located under the bench and secure it with the supplied hose clamp.*

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*ALL MAINTENANCE MUST BE PERFORMED WITH THE UNIT OFF AND THE POWER PLUG REMOVED FROM THE OUTLET!*

*The hood is an easy-to-use appliance requiring little maintenance. Following are a few directions:*

*The appliance filter located below the exhauster must be periodically cleaned, depending on the working hours, but at least every 500 working hours.*

*To carry out the filter cleaning, undo the hooks on the sides, extract the filter holder and retainer.*

*Remove the filter and clean it.*

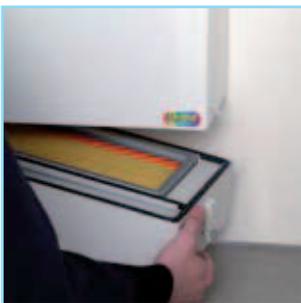
*CAUTION: while reassembling, do not forget to reinsert the filter retainer.*

*In any case, the filter must be replaced after 1 year of use.*

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*To clean or to replace the filter, undo the hooks on the sides of the hood.*

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*Grasp and extract the filter holder as shown in the picture.*

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*Remove the metal filter retainer. CAUTION! Do not forget to reinsert it after cleaning or replacing the filter.*

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*Remove the filter as shown in the picture and clean or replace it. To reassemble, repeat the previous steps in reverse order.*

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**IRIDE INTERNATIONAL S.R.L.**

## DECLARATION OF CONFORMITY

We, IRIDE INTERNATIONAL S.r.l. – furniture for dentist offices and dental laboratories, with registered headquarters in Via San Giuseppe 5 - 33082 Azzano Decimo - Località Tiezzo - PN:

**HEREBY DECLARE UNDER OUR OWN RESPONSIBILITY THAT THE PRODUCT “BENCH FUME HOOD” TO WHICH THIS DECLARATION RELATES WERE DESIGNED AND BUILT IN EXCLUSIVE MANUFACTURING IN CONFORMITY WITH THE FOLLOWING HARMONISED STANDARDS:**

**IEC EN 60601-1(IEC 65-5) of 1991 and its 1994 amendments**

**IEC EN 60204-1(IEC 44/5 and 44/6)**

**AND THE REQUIREMENTS OF THE DIRECTIVE:**

**2006/42/CE**

**2006/95/CE**

**2004/108/CE**

**IRIDE INTERNATIONAL S.R.L.**

**IRIDE INTERNATIONAL s.r.l.**

Via S. Giuseppe 5 | 33082 Azzano Decimo - PN

tel. +39 0434 633242 | fax +39 0434 632546

welcome@irideinternational.com | www.irideinternational.com